



Rapid Characterisation of Battery Materials using Raman Spectroscopy

Tarun Sharda, M.Sivakumar, N.Sathish

TechnoS Instruments, Jaipur, India.

CSIR-AMPRI, Bhopal, India.

INTRODUCTION

Greatest challenge of current era is to meet the energy demand by providing the proper energy supply which is hampered by increasing oil shortage and rising pollution level due to excessive coal usage. Relying on the batteries is one of the best solutions to address the existing problems and to meet the growing energy demand. Materials plays a vital role in deciding the efficiency and lifespan of the manufactured batteries. Fundamental understanding of materials and its characterization is very important in taking the technology forward in a quick phase. Though different techniques are available for the identification and characterization of battery materials, Raman spectroscopy is one of the quickest techniques among them. Also, it is essential to extract the data from the working electrodes in-situ to understand the interactions of materials which is possible by Raman Spectroscopy. The objective of the current study is the quick characterization of the battery materials used as electrodes, electrocatalysts and identification of number of layers of reduced graphene oxide (rGO) layers by mapping using Raman spectroscopy.

MATERIAL AND METHODS

Raman Spectroscopy of the battery materials was performed using an IndiRAMTMCTR-500 Raman spectrometer, TechnoS Instruments Ltd., Jaipur, India. A laser excitation source of 532 nm emission wavelengths was used. Different industrial grade graphene samples, anode and electrocatalyst materials were characterized.



IndiRAM CTR-500 Raman spectrometer

RESULT AND DISCUSSION

Figure 1(a) shows the Raman spectra acquired from different industrial grade Graphene samples showing G and D band. Figure 1(b) shows the spectra of ZnTiO_3 , $\text{Sr}_2\text{TiMnO}_6$, and $\text{CaCu}_3\text{Ti}_4\text{O}_{12}$ respectively. In ZnTiO_3 spectrum, 311 and 446 cm^{-1} are due to the oxygen vibration of Zn–O–Zn, 610 cm^{-1} corresponds to vibrations of Ti–O–Ti[1]. Convoluted asymmetric peak around 500 cm^{-1} (475 and 523 cm^{-1}) in $\text{Sr}_2\text{TiMnO}_6$ corresponds to the T_{2g} vibration modes [2]. The Raman lines of $\text{CaCu}_3\text{Ti}_4\text{O}_{12}$ at 446, 510 and 572 cm^{-1} correspond to Ag symmetry and Fg symmetry respectively [3]. Fig 1 (c) shows mapping of reduced GO in which layer numbers are directly related to the IG ratios and the corresponding intensities and peak positions are tabulated in

fig1 (a)

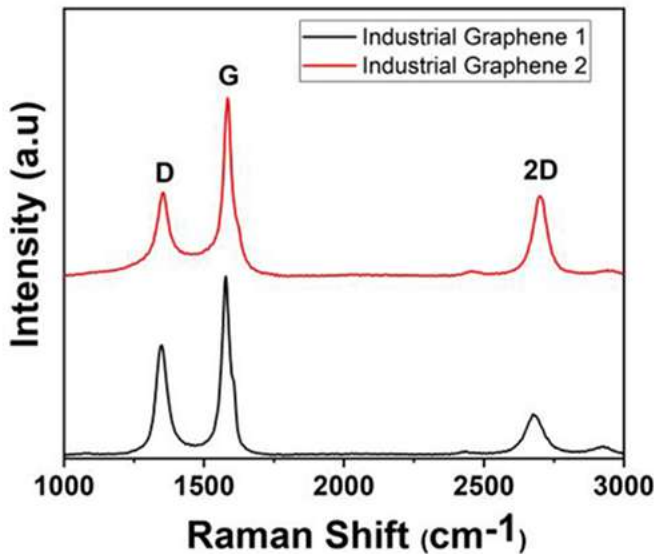


fig1 (b)

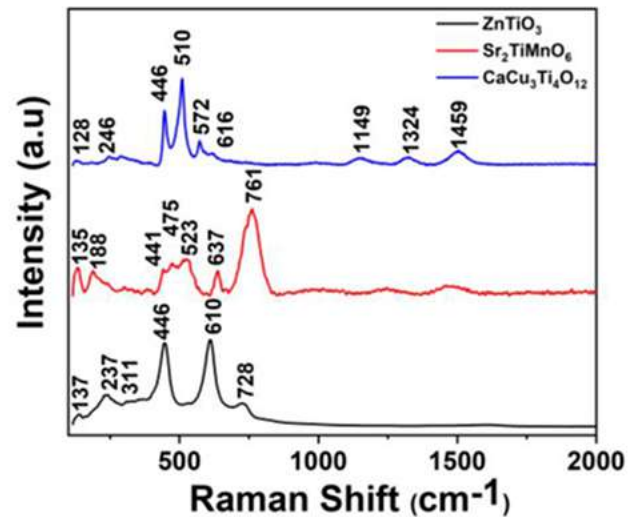


fig1 (c)

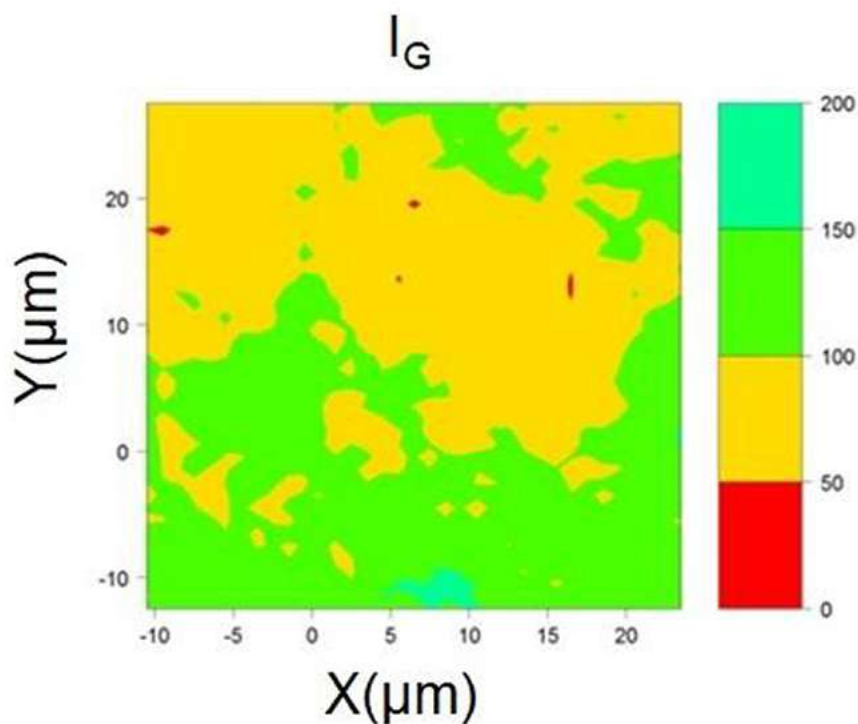


Table:1 Intensity and Peak positions

Region	I_D/I_G	I_G	G	D	I_G Ratio
Yellow	1	4816	1580	1348	2
Green	1	6978	1579	1346	2.9

CONCLUSION

In this application note, quick characterization of battery materials has been carried out by Raman Spectroscopy.

REFERENCES

1. Arin, J., Thongtem, S., Phuruangrat, A. & Thongtem, T. Characterization of ZnO–TiO₂ and zinc titanate nanoparticles synthesized by hydrothermal process. Research on Chemical Intermediates 43, 3183–3195 (2017).
2. Murugesan, G., Nithya, R. & Kalainathan, S. Colossal dielectric behaviour of Sr₂TiMnO₆- δ single crystals. J Cryst Growth 530, (2020).
3. Kolev, N. et al. Raman spectroscopy of (formula presented). Phys Rev B Condens Matter Mater Phys 66, 1–4 (2002).

OUR PRODUCTS



IndiRAM™ CTR Series



IndiRAM™ CTR for
Quantum Characterizer



IndiRAM™ CTR Series
Chiral Raman Spectrometer



IndiRAM™ Edu



IndiRAM™ CTR
P-Series



IndiRAM™ Portable
System



Diamond Detection
System



IndiRAM™
Handy Series



IndiRAM™ CTR-
Mini Series

